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Vivapore 2, 5, 10 | 20 ml C€

Directions for Use

Intended only for in-vitro diagnosis

Vivapore Concentrators offer exceptional ease of use and flexibility for the concentration and/or purification of macromolecules from dilute solutions. Depending on configuration and reservoir volume of the Vivapore selected, starting volumes may vary from 0.5 to 20 ml and up to 750 concentration may be achieved.


Operation

1. Concentrating macromolecules

a. Insert Vivapore 5 or 10 | 20 into stand. Vivapore 2 is self-supporting.

b. Pipette sample through aperture at the top of the device. Device can be left unattended until desired concentration is achieved. Solvent and micromolecules are pulled through the membrane by a high-capacity absorbent; a dead stop feature prevents the sample from concentrating to dryness.

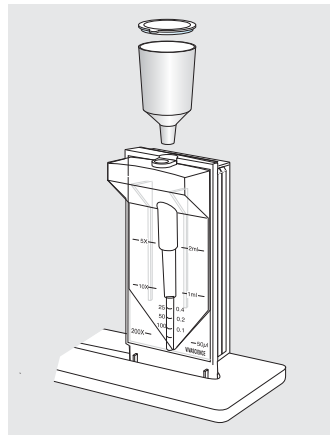c. Once the desired volume is achieved, the concentrate is withdrawn using a Pasteur, thin plastic or gel loader type pipette. The sample is now ready for further analysis.

2. 20 ml capacity expansion for the VP10 | 20

Sample capacity of the VP10 | 20 can be increased to 20 ml by use of the ancillary reservoir. First pipette 10 ml into the Vivapore, then lock the ancillary reservoir to the top opening and add up to an additional 10 ml of sample.

3. Desalting with either Vivapore 2 or Vivapore 10 | 20


Reservoir expansion or desalting by constant volume diafiltration. Both the VP10 | 20 and Vivapore2 offer a unique ancillary reservoir feature that allows reservoir expansion or one-step purification and concentration of protein samples.

In this procedure, new buffer solution replaces the original solvent at a rate equal to the speed of filtration. With a fresh buffer solution volume that is five times greater than the sample in the concentrator, 99% of salts will be removed.

A. Pipette up to 2 ml of sample to be purified through aperture at the top of the device.

B. Firmly seat a 10 ml plastic pipette or the ancillary reservoir on the conical entry port of the concentrator. Fill with up to 5 times the volume in the concentrator using deionized water or fresh buffer solution. Device can be left unattended whilst the new buffer replaces the solvent in the original solution. Once buffer in the pipette is exhausted, the remaining solution will continue to concentrate to the level desired.

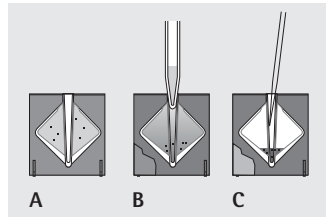C. Collect purified and concentrated sample with a pasteur or long gel loader type pipette.

Operating Tips

Improving speed of concentration

Speed of filtration is affected by several parameters including temperature, pH and protein concentration. Whilst Vivapore concentrators will provide rapid filtration in most environments, the following suggestions may be helpful in optimising speed.

Filtration speed will increase proportionally to ambient temperature. Should you require faster concentration, place the concentrator near a source of heat.

An acid sample with a pH of less than 5 will take longer to concentrate than a neutral sample. Adjustment to a physiological pH will result in faster filtration.

Suspended particles will tend to foul the filter element and slow filtration speed. Prefiltration with a syringe filter will clarify the sample and result in faster filtration speed and improved analytical results following concentration.

Initial protein concentration levels will have a significant effect on concentration speed. Whilst a highly dilute sample will concentrate rapidly, once macromolecule concentration exceeds 2%, the speed of filtration will rapidly decrease.

Concentrations above 5% are not practical with Vivapore concentrators. (For high levels of macromolecular concentration, use Vivaspin centrifugal concentrators.)

Recovering dry samples

Vivapore concentrators have an impermeable concentrate pocket (dead stop) which impedes concentration to dryness. However, if the concentrate inadvertently remains too long in the concentrator, the remaining solvent will eventually evaporate and the sample may go to dryness. Should this occur, proteins may be returned to solution by pipetting approximately 100 µl of buffer in and out of the concentrate pocket several times.

Storage

Vivapore devices should be stored in a dry environment at 4 to 30°C.

Disposal

Used units may be incinerated or autoclaved, the stands are reusable for further concentrations.

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* with additional reservoir

Table 2: Performance characteristics

Product	Time to concentrate up to 10x [min.]				Concentrate recovery %			
	VP2	VP5	VP10 20	VP10 20*	VP2	VP5	VP10 20	VP10 20*
Start volume	2 ml	5 ml	10 ml	20 ml	2 ml	5 ml	10 ml	20 ml
Cytochrome c (12,600 MW) 7,500 MWCO PES	0.25 mg/ml 35	0.25 mg/ml 35	0.25 mg/ml 75	0.1 mg/ml 150	0.25 mg/ml 90 %	0.25 mg/ml 90 %	0.25 mg/ml 90 %	0.1 mg/ml 92 %
BSA (66,000 MW) 7,500 MWCO PES	25	30	55	115	90 %	92 %	92 %	92 %
IgG (160,000 MW) 7,500 MWCO PES	35	40	70	160	76 %	75 %	77 %	78 %
	Time to concentrate up to 50x [min.]				Concentrate recovery %			
Cytochrome c (12,600 MW) 7,500 MWCO PES	65	70	160	-	91 %	88 %	90 %	-
BSA (66,000 MW) 7,500 MWCO PES	45	50	105	218	90 %	90 %	92 %	94 %
IgG (160,000 MW) 7,500 MWCO PES	50	65	140	290	53 %	65 %	74 %	70 %

* with additional reservoir

Table 3: Ordering information

	Pack size	Prod. no.	
Vivapore 2 Expandable to 15 ml with pipette reservoir 7,500 MWCO PES	30	VP0201	
Vivapore 5 Includes stand and recovery pipettes 7,500 MWCO PES	4	VP0503	
7,500 MWCO PES	30	VP0501	
Requires stand 7,500 MWCO PES	100	VP0502	
Vivapore 10 20 Includes stand and recovery pipettes 7,500 MWCO PES	4	VP2003	
7,500 MWCO PES	30	VP2001	
Requires stand 7,500 MWCO PES	100	VP2002	
Vivapore accessories			
Disposable stands for 4 units	6	VPA002	
Pipette reservoir (Vivapore 2)	50	VPA004	
Plastic recovery pipettes (Vivapore 10 20)		100	VPA005
10 ml expansion reservoir (Vivapore 10 20)	10	VPA006	
Plastic recovery pipettes (Vivapore 5)	100	VPA007	

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany
Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

The information in this publication is believed to be accurate, and is offered in good faith for the benefit of the user. As we cannot anticipate or control conditions of product application, we do not warrant suitability or favourable results, nor that use will not infringe on any patents. Products are furnished upon the condition that the user assumes all risks and liabilities and that neither seller or manufacturer shall be liable for any loss or damage, direct or sequential arising from the use or inability to use the product.

Specifications subject to change without notice. Printed and copyrighted by Sartorius · W405.21
Publication No.: SLU6098-p08023